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3.6 Hamilton Syringes 3. Accessories, Spare Parts, Syringes etc.

Choosing the Right Syringe

- **Which termination works best for your application?**
- **Luer Tip Cemented** - ground glass tip with a cemented needle ; not recommended for use with halogenated solvents, such as MeCl₂
- **Teflon® Luer Lock** - can be used with luer (Metal or Kel-F®) hub needles, Teflon® tubing assemblies and connectors can be used, offer flexibility
- **Cemented** - not recommended for use with halogenated solvents, such as MeCl₂
- **Luer Tip** - ground glass tip can be used with most hypodermic needles; Kel-F® hub needles, Teflon® tubing assemblies and connectors can also be used; offers flexibility
- **Removable** - advantageous for laboratories with multiple users; economical since you can replace the needle without having to purchase a new syringe

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3.6 Hamilton Syringes

3. Accessories, Spare Parts, Syringes etc.

Selecting the Right Syringe

Tips to Assure Accuracy and Precision

- When using a Hamilton syringe, grasp only the syringe flange and plunger button. By doing so, variations in liquid measurement due to body heat are avoided.
- Pump the plunger with the syringe needle immersed in the fluid to be transferred. This will expel any trapped air in the needle and syringe. Minimize the use of a dry syringe.
- If the plunger is accidentally removed from the syringe barrel wipe it carefully with a lint-free tissue. Reinsert the plunger into the barrel and pump deionized water or acetone through the needle and syringe. In the case of GASTIGHT® plunger, re-wet the Teflon® PTFE plunger tip prior to reinserting it into the barrel. Avoid touching the plunger since any abrasions, scratches, or oil from one's fingers will often interfere with the proper plunger operation.

Syringe Lubrication

A clean syringe does not require any lubrication grease. Grease may cause a variety of problems including: sample cross contamination, seizing of the plunger in the barrel or barrel damage.

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Choosing the Right Needle Style

Point Style 3

blunt needle point for use with HPLC injection valves and for sample pipetting.

Point Style AS

special conical style needle point designed to withstand the demands of multiple injections; exclusively used on auto-sampler syringes.

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Several different needle point styles are offered on Hamilton syringe and needle products depending upon the application. Syringes and needles manufactured by Hamilton Company are intended for scientific research and laboratory use only and are not intended for human in vivo use.

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HAMILTON
the measure of excellence

3.6 Hamilton Syringes

3. Accessories, Spare Parts, Syringes etc.

Explanation and Abbreviation

Syringe Terminations

N, NR (Cemented Needle, Cemented Needle Rheodyne)

Needles are cemented into the glass syringe barrel at a point corresponding to the zero graduation mark. NR stands for syringes with Rheodyne specified needles.

LT (Luer Tip)

Ground glass syringe barrel with a male luer taper accepts most hypodermic needles. Use Kel-F® hub needles and connectors for a tight seal.

RN, RNR (Removable Needle, Removable Needle Rheodyne)

Removable needles seat precisely to the zero graduation mark of the syringe. Allows the use of different specification needles on the same syringe barrel. RNR stands for syringes with Rheodyne specified removable needles.

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3.6 Hamilton Syringes

3. Accessories, Spare Parts, Syringes etc.

...for Rheodyne, Valco™, Beckman®/Altex, SSI, KNAUER Injection Valves

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3.6 Hamilton Syringes

3. Accessories, Spare Parts, Syringes etc.

Syringe Cleaning Tips

Cleaning Tips

- The life of your MICROLITER™, MODIFIED MICROLITER™ or GASTIGHT® syringe is directly related to its cleanliness!
- To clean syringes, it is best to use solvents know to be effective in solvating the sample and preferably are non-alkaline, non-phosphate and non-detergent based. A biodegradable, non-phosphate, organic Cleaning Concentrate is available from Hamilton; order part number HA 18311.
- High quality water and acetone prove to be good rinses.
- To clean the plunger, remove it from the syringe barrel and gently wipe with a lint-free tissue. Reinsert the plunger into the barrel and pump deionised water or acetone through the needle and syringe. Air dry the syringe for storage.

Note: When reinserting a Teflon® -tipped plunger into a syringe barrel, wet (lubri cate) the tip with deionised water or another solvent compatible with the sample.

Sterilizing Syringes

Sterilize your Hamilton syringe using an appropriate chemical sterilizer. Rinse the syringe thoroughly after sterilization. In general, Hamilton does not recommend autoclaving syringes. Stainless steel expands faster than glass upon heating. Heating a syringe with a cemented, stainless steel needle may cause the glass barrel to split. Heating a syringe with cemented stainless steel hubs may result in adhesive deterioration and leakage. It is the responsibility of the user to determine the suitability of application and material compatibility of the products based on published specifications.

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Needle Cleaning Kit

- Contains a selection of various diameter tungsten wires for cleaning plugged needles
- Biodegradable cleaning concentrate helps to remove tenacious residues

Bestell Informationen / Ordering Information

Part Number	Description																					
HA 76620	Needle Cleaning Kit (includes cleaning wires and 70 ml of cleaning concentrate listed below)																					
	<table border="1"> <thead> <tr> <th>Cleaning Wire Part Number</th> <th>Cleaning Wires (178 mm in length)</th> <th>Use to Clean Needle Gauges</th> </tr> </thead> <tbody> <tr> <td>HA 18306</td> <td>0.076 mm O.D., 10/pk</td> <td>23s, 26s, 31-33</td> </tr> <tr> <td>HA 18300</td> <td>0.089 mm O.D., 10/pk</td> <td>26s, 31-33</td> </tr> <tr> <td>HA 18301</td> <td>0.126 mm O.D., 10/pk</td> <td>22s, 25s, 28-30</td> </tr> <tr> <td>HA 18302</td> <td>0.167 mm O.D., 10/pk</td> <td>27</td> </tr> <tr> <td>HA 18303</td> <td>0.207 mm O.D., 10/pk</td> <td>24-26</td> </tr> <tr> <td>HA 18304</td> <td>0.306 mm O.D., 10/pk</td> <td>22, 23 (and larger)</td> </tr> </tbody> </table>	Cleaning Wire Part Number	Cleaning Wires (178 mm in length)	Use to Clean Needle Gauges	HA 18306	0.076 mm O.D., 10/pk	23s, 26s, 31-33	HA 18300	0.089 mm O.D., 10/pk	26s, 31-33	HA 18301	0.126 mm O.D., 10/pk	22s, 25s, 28-30	HA 18302	0.167 mm O.D., 10/pk	27	HA 18303	0.207 mm O.D., 10/pk	24-26	HA 18304	0.306 mm O.D., 10/pk	22, 23 (and larger)
Cleaning Wire Part Number	Cleaning Wires (178 mm in length)	Use to Clean Needle Gauges																				
HA 18306	0.076 mm O.D., 10/pk	23s, 26s, 31-33																				
HA 18300	0.089 mm O.D., 10/pk	26s, 31-33																				
HA 18301	0.126 mm O.D., 10/pk	22s, 25s, 28-30																				
HA 18302	0.167 mm O.D., 10/pk	27																				
HA 18303	0.207 mm O.D., 10/pk	24-26																				
HA 18304	0.306 mm O.D., 10/pk	22, 23 (and larger)																				
HA 18311	Cleaning Concentrate, 500 ml (not included in cleaning kit)																					

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Cleaning Needles

Should a needle become clogged, do not attempt to clean it by forcing liquid or compressed air through the syringe. Excessive pressure will split the glass barrel. Order the Needle Cleaning Kit, part number HA 76620, to prevent clogging of needles. Remove the plunger from the syringe before cleaning the needle. Cleaning wires are used to dislodge any foreign material or residue build-up. Use deionised water or acetone to thoroughly rinse the syringe.

Syringes

Notizen

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Notes

3. Accessories, Spare Parts,
Syringes etc.

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