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The Headspace instruments on the market require a broad variety of vials and closures. To facilitate the identification of the right vial, indications on the instrument suitability are given at the bottom of each table.

The right choice of closure might depend on the instrument or application dependent. Therefore please study the individual features of the various types of caps in chapter 5.9.3 as well as the properties of the liners at the top of each seal chapter. In case of questions, you are always welcome to contact us for advice.

A Screw Neck Vial for PERKIN ELMER with PP Screw Seals as well as Screw Neck Vials ND18 with Magnetic Screw Seals for CTC are novelties in the crimp neck dominated Headspace market. However, especially for the hard to crimp magnetic caps, screw closures are a convenient alternative.

IMPORTANT SAFETY TIP:

As Headspace vials have to withstand internal pressure, almost all vials on the next page have a wall thickness of 1.2 mm. Thus it is guaranteed that the vial will not burst. Contrary to a widely spread opinion the seal and not the vial represents the weakest part of the whole system. Under pressure the septa will bulge against the Aluminium cap with such a force that the cap is torn apart. By testing is verified that e.g. the Perkin Elmer vial LP20 09 0297 can easily withstand 10 bars or more while the seal is torn apart at around 10 bars when no pressure release system - like our Headspace Cap - is used.

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5. HPLC Vials

Headspace Vials might differ in:

- Volume (5 ml, 10 ml, 20 ml)
- Rounded or flat bottom
A rounded bottom is more sturdy and thus more resistant to the high pressure within the vial during the heating process. Furthermore it slides more easily into the heating block when being transported by a magnet.
A flat bottom might be necessary when vials have to run within the instrument on a slightly downwards tendency.
- Bevelled top or flat DIN Crimp Neck / Screw Neck
Bevelled top might be required for some special closure systems (PERKIN ELMER), however a liner has more surface to lie on with a flat DIN Crimp Neck. The more surface for the liner to rest on, the tighter the seal.
- Length of the Neck (instrument specific)
- Clear / Amber Glass
- With / without label and filling lines
- Upon request with a barcode label

Headspace Closures might differ in:

- Type of Cap
Centre Hole Crimp Cap, Headspace Cap, Centre Tear Off Cap, Complete Tear Off Cap, Magnetic Cap, SPME Cap, Magnetic Screw Cap, PP Screw Cap, PE Push-on Cap.
Screw Caps are a novelty in the crimp neck dominated Headspace area. However, they represent a ready to use, convenient solution, that does not require any additional tools (crimpers, decappers). Thus samples can be taken and sealed out in the field without the necessity of sample transfer later in the lab. The magnetic Screw Seals can be used universally for Headspace as well as for SPME.
- Butyl, Butyl/PTFE, Pharma-Fix-Liner (Butyl/PTFE), Silicone/PTFE, Silicone/Aluminium Foil, Viton, Natural Rubber/TEF.
Besides the material the liners might differ in thickness, hardness (° shore A), colour, type of PTFE lamination, grade of Silicone (UltraClean™)

5.9 Headspace ND20 (ND18)

5. HPLC Flaschen

5.9.1 Headspace-Flaschen ND18 und ND20

Headspace-Flaschen ND18 und ND20

LP20 09 0342	LP20 09 0801	LP20 09 1405	LP20 09 0802	LP20 09 0795	LP20 09 0297 LP20 09 1223	LP20 09 0440
5.0 ml Headspace Vial 38.2 x 22 mm, clear glass, 1st hydrol. class, rounded bottom	5.0 ml Crimp Neck Vial 38 x 20 mm, clear glass, 1st hydrol. class, flat bottom	10 ml Headspace Vial 46 x 22.5 mm, clear glass, 1st hydrol. class, DIN -Crimp Neck, rounded bottom	10 ml Crimp Neck Vial 54.5 x 20 mm, clear glass, 1st hydrol. class, flat bottom	10 ml Headspace Vial 46 x 22.5 mm, clear glass, 1st hydrol. class, DIN-Crimp Neck, long neck, flat bottom	20 ml Headspace Vial 75.5 x 23 mm, clear/amber glass, 1st hydrol. class, rounded bottom,	20 ml Headspace Vial 75.5 x 23 mm, clear glass, 1st hydrol. class, rounded bottom, with label + filling lines
100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box
Perkin Elmer	Varian	Carlo Erba, CTC, Fisons, Varian (CP)	Varian	Carlo Erba, Dani, Fisons, Agilent	Perkin Elmer	Perkin Elmer

HPLC Headspace ND20 (ND18)

5.9 Headspace ND20 (ND18)

5. HPLC Vials

5.9.1 Headspace Vials ND18 and ND20

Headspace Vials ND18 and ND20

HPLC Headspace ND20 (ND18)

5.9 Headspace ND20 (ND18)

5. HPLC Flaschen

5.9 Headspace ND20 (ND18)

5. HPLC Vials

5.9.1 Headspace-Flaschen ND18 und ND20

5.9.1 Headspace Vials ND18 and ND20

Headspace-Flaschen ND18 und ND20

Headspace Vials ND18 and ND20

LP20 09 0286	LP20 09 0796	LP20 09 0873	LP20 09 1222	LP18 09 0864	LP18 09 1306 LP18 09 1310	LP18 09 1307 LP18 09 1311
20 ml Crimp Neck Vial 75.5 x 23 mm, clear glass, 3rd hydrol. class, flat bottom	20 ml Headspace Vial 75.5 x 22.5 mm, clear glass, 1st hydrol. class, DIN-Crimp Neck, long neck, rounded bottom	20 ml Headspace Vial 75.5 x 22.5 mm, clear glass, 1st hydrol. class, DIN-Crimp Neck, long neck, rounded bottom	20 ml SPME Vial 75.5 x 22.5 mm, clear glass, 1st hydrol. class, rounded bottom, special crimp neck	20 ml Headspace Vial 75.5 x 23 mm, clear glass, 1st hydrol. class, rounded bottom, with screw thread ND18	10 ml Screw Neck Vial ND18, 46 x 22.5 mm, clear/amber glass, 1st hydrol. class, rounded bottom	20 ml Screw Neck Vial ND18, 75.5 x 22.5 mm, clear/amber glass 1st hydrol. class, rounded bottom
100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box	100 pcs./PP-Box
Perkin Elmer	Carlo Erba, Dani, Fisons, Agilent	CTC PAL (Varian, Gerstel, Atas, Shimadzu)	SPME Vial for CTC PAL	Perkin Elmer	CTC PAL (Varian, Gerstel, Atas, Shimadzu)	CTC PAL (Varian, Gerstel, Atas, Shimadzu)

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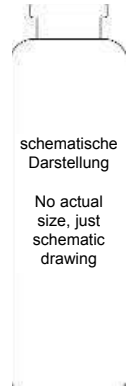
Sonstige Rollrandflaschen ND20

LP20 09 0289

50 ml
Crimp Neck Vial,
101 x 31 mm,
clear glass
1st hydrol. class

100 pcs./PP-Box

5.9 Headspace ND20 (ND18)

5. HPLC Vials

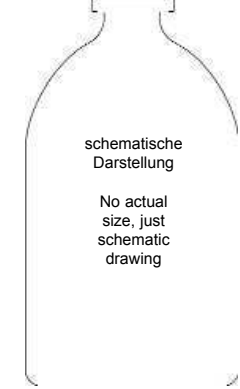
5.9.2 Other Crimp Neck Vials ND20

Other Crimp Neck Vials ND20

LP20 09 0343

100 ml
Crimp Neck Vial,
95 x 52 mm,
clear glass
3rd hydrol. class

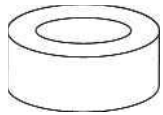
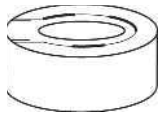
72 pcs. shrink-wrapped

Any 20 ml Headspace-Vial can be obtained with the appropriate seal as a 2in1 KIT!

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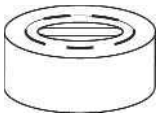
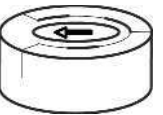
5.9.3 Aluminium Crimp Seals ND20

A broad selection of different liners can be obtained in the following caps:

	Mittelabrisskappe <i>Centre Tear-Off Cap</i>	Ganzabrisskappe <i>Complete Tear-Off Cap</i>
	1:1 	1:1 
	Clear lacquered, red, blue, gold, green	Clear lacquered, red, blue, gold, green

Any type of cap can ONLY be obtained in combination with a liner or a stopper!

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5.9.3 Aluminium Crimp Seals ND20

5.9.3.1 Butyl Seals

- Temperature resistant from -40 °C up to 120 °C
- Due to the missing PTFE lamination only suitable for uncritical analysis
- Low-cost product

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5.9.3.2 Butyl/PTFE Seals

- Temperature resistant from -40 °C up to 120 °C
- Completely laminated with PTFE

Packaging Unit: 100 pcs./PE-Bag